



Echelon Biosciences Inc.
675 Arapeen Drive, Suite 302
Salt Lake City, UT 84108
Telephone 866-588-0455
Fax 801-588-0497
echelon@echelon-inc.com
www.echelon-inc.com

Technical Data Sheet

For research use only
Not intended or approved for
diagnostic or therapeutic use.

Sphingo Strips™

Product Number: S-6000

Product Description:

Sphingo Strips™ are 2 x 6 cm hydrophobic membranes that have been spotted with 15 different lipids at 100 pmol per spot. These membranes can be used to determine lipid-protein interactions, through a simple protein-lipid overlay experiment. This allows researchers a convenient way to determine if their protein of interest interacts with one or more of the bound lipids.

Storage:

Store at 2-8 °C. Product is moisture and light sensitive.

Format:

The membrane has a diagonal cut on its top left corner and is spotted with Xylene Cyanol FF (blue) in the bottom right blank corner to assist in orientation of the strip. Ponceau S staining (pink) was added to the lipid spots. See template below for location of lipids. All of the lipids are long chain (> diC16) highly pure synthetic analogs. For more information, on the lipids spotted on the membrane, please visit our website and search the catalog numbers provided in the figure below.

Membrane Template:

Sphingosine (cat # S-1000)	○ ○	Monosialoganglioside-GM1
Sphingosine-1-phosphate (S1P, cat # S-2000)	○ ○	Disialoganglioside-GD3
Phytosphingosine	○ ○	Sulfatide
Ceramide	○ ○	Psychosine
Sphingomyelin (SM)	○ ○	Cholesterol (cat # L-6012)
Sphingosylphosphorylcholine (SPC)	○ ○	Lysophosphocholine (LPC, cat # L-1518)
Lysophosphatidic Acid (LPA, cat # L-0200)	○ ○	Phosphatidylcholine (PC, cat # L-1116)
Myriocin	○ ●	Blue Blank

*Final concentration of 0.1% (v/v) Ponceau S was added for accuracy during spotting.

Suggested Usage:

Visit our website www.echelon-inc.com. At the bottom of the each product's webpage is our general Protocol "Protocol_Strip_Array" for use with product numbers: P-6001, P-6100, P-6002, P-6003, S-6000, and S-6001. Also please refer to our FAQ "Frequently Asked Questions" document.

References:

1. Dowler, S., Currie, R.A., Downes, P.C., and Alessi, D.R. DAPP1: a dual adaptor for phosphotyrosine and 3-phosphoinositides *Biochemical Society J.* 342, 7-12 (1999).
2. Dowler, S., Kular, G., and Alessi, R.D., Protein lipid overlay assay, *Sci STKE*, 2002, L6. (2002).
3. Rodriguez-Asiain A, Ruiz-Babot G, Romero W, Cubi R, Erazo T, Biondi RM, Bavascas JR, Aquilera J, Gomez N, Gil C, Claro E, Lizcano JM. Brain specific kinase-1 BRSK1/SAD-B associates with lipid rafts: modulation of kinase activity by lipid environment. *Biochem Biophys.* 2011 Dec.
4. Sawa CG, Fernandes da Costa SP, Bokori-Brown M, Naylor CE, Cole AR, Moss DS, Titball RW, Basak AK. Molecular architecture and functional analysis of NetB, a pore-forming toxin from *Clostridium perfringens*. *J Biol Chem.* 2013 Feb.

Echelon Biosciences products are sold for research and development purposes only and are not for diagnostic use or to be incorporated into products for resale without written permission from Echelon Biosciences. Materials in this publication, as well as applications and methods and use, may be covered by one or more U.S. or foreign patents or patents pending. We welcome inquiries about licensing the use of our trademarks and technologies at busdev@echelon-inc.com.