


# Mobicol "F" with fixed outlet plug, inserted 35 µm filter and 2 different screw caps

## Product Information Sheet # M105035F

### SUMMARY

shipped at RT; store at RT

For research use only

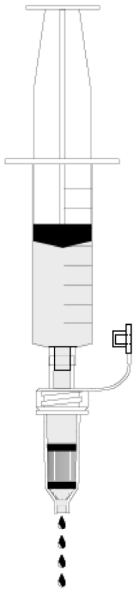

Syringe connected to Luer-lock cap


Mobicol "F" is a 1 ml column with particular handling advantages.

Mobicol "F"

- is compatible to the laboratory standard,
- can be centrifuged in minifuges,
- has a Luer-lock adapter,
- smallest and largest volumes can be treated easily.



Upper filters of the pore sizes 10 µm, 35 µm or 90 µm are available for Mobicols. The upper large filter fits tightly over the whole length of the column body.

Mobicols can be closed tightly with a screw-on cap and the reversed side of the snap-off plug at the outlet. The volume is 700 µl. Mobicols can be autoclaved at 110 °C for 10 minutes.

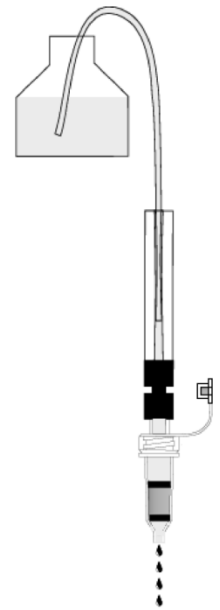

Adapter M3001: tubing to Mobicol

### Mobicol "F"

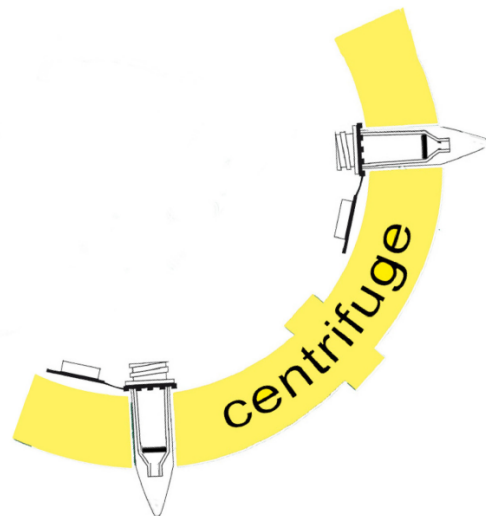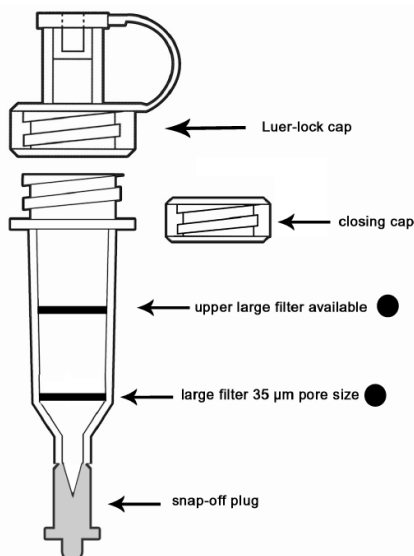

Mobicols can be centrifuged (fit into 1.5 ml or 2 ml tubes)

# Mobicol "F" with fixed outlet plug, inserted 35 µm filter and 2 different screw caps

## Product Information Sheet # M105035F



The column M105035F is supplied with two screw caps (Luer-lock cap and screw cap) and a snap-off plug for the outlet. Upper filters with different pore sizes (10 µm, 35 µm and 90 µm) are separately available, so you can use the filter of your choice. With this features the Mobicol "F" can be used for a variety of applications.

### Order Information, Shipping and Storage

Order#	Product	Quantity
M105035F	Mobicol "F" with fixed outlet plug, inserted 35 µm filter and 2 different screw caps	50
M1050	Mobicol "F" with fixed outlet plug and 2 different screw caps, without filters	50
M1053	Mobicol "F" with fixed outlet plug and screw cap, without filters	50
M2210	Filter (large) 10 µm pore size, for Mobicol M1002, M1003, M1050 & M1053	50
M523515	Filter (large) 35 µm pore size, for Mobicol M1002, M1003, M1050 & M1053	50
M2290	Filter (large) 90 µm pore size, for Mobicol M1002, M1003, M1050 & M1053	50
shipped at RT; store at RT		

### Column Accessories

Order#	Product	Quantity
M3001	Tubing-to-Luer-lock adapter (material outside: metal; inside: glass and PTFE; with 1 m PTFE tubing) for Mobicols or laboratory columns	1
M3002	Luer adapter, (M+F) 10 each	20
M3003	Luer adapter for the top-cap (M)	20
M3004	Luer adapter for outlet (F)	20
M3006	Stopcocks, one-way, Luer-lock	4
M3007	Luer-lock adapter (M/M)	1
shipped at RT; store at RT		

### Contact and Support

MoBiTec GmbH ● Lotzestrasse 22a ● D-37083 Goettingen ● Germany

#### Customer Service – General inquiries & orders

phone: +49 (0)551 707 22 0  
fax: +49 (0)551 707 22 22  
e-mail: [order@mobitec.com](mailto:order@mobitec.com)

#### Technical Service – Product information

phone: +49 (0)551 707 22 70  
fax: +49 (0)551 707 22 77  
e-mail: [info@mobitec.com](mailto:info@mobitec.com)

**MoBiTec in your area:** Find your local distributor at [www.mobitec.com](http://www.mobitec.com)