

Dimensions of the Adapter*12,75 x 8,55 x 14 mm*


	Position 1	Position 2	Position 3	Position 4
X	10 mm	37,25 mm	64,5 mm	91,75 mm
Y	4 mm	4 mm	4 mm	4 mm

The coordinates of the different patterns can be found in the file: "Specifications of the DMA". (Download at: <https://aquarray.com/product-information>) The coordinates given are for position 1. Please adjust for the other positions by adding the values for x in the above table.


Loading of the DMA slide into the Adapter

To address the full field of the pattern, the adapter has to be used upside down.


Place the adapter on the table with position 1 on the left side.

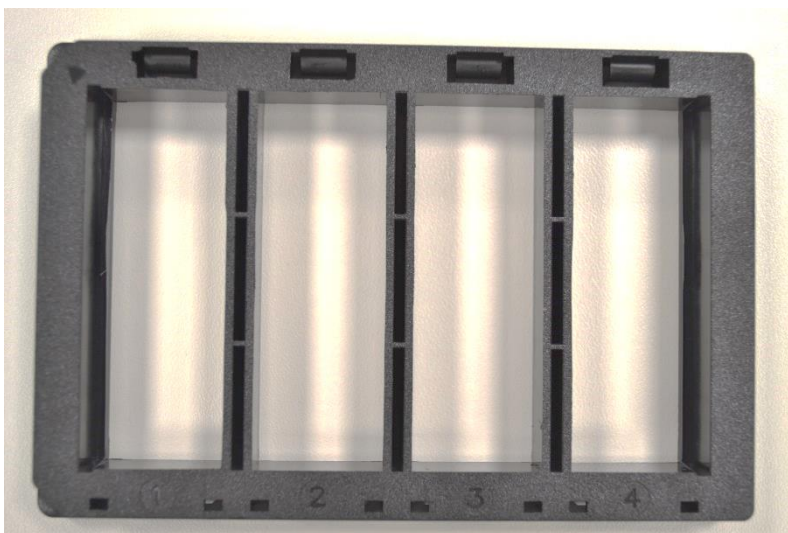

To load the DMA Slide in the adapter, turn the adapter upside down. Position 1 should stay on the left side.

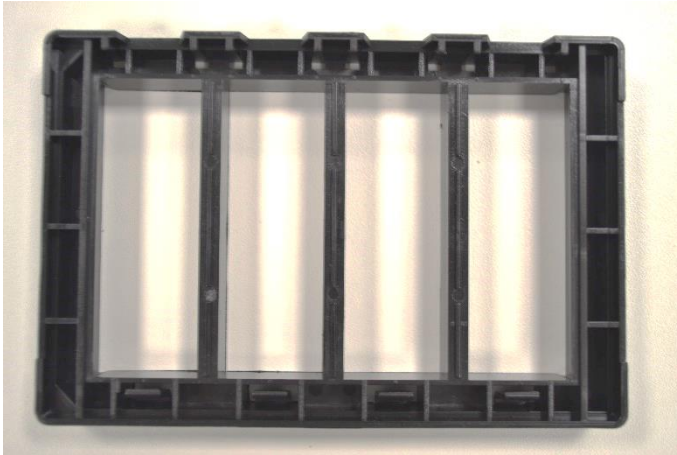

Place a DMA in position 1 elow the two holding pins. Push the DMA slide to the left side using tweezers. Carefully push from the sides to prevent scratches on the surface.

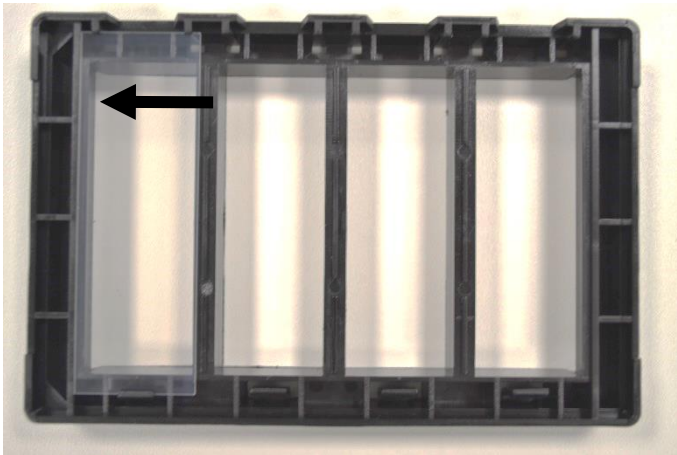

Exception:

For DMA slides with the product number *-102* and *-401* please push the slide to the right side using tweezers.

