ProFoldin



10 Technology Drive, Suite 40, Number 188 Hudson, MA 01749-2791 USA Tel: (508) 735-2539 FAX: (508) 845-9

Tel: (508) 735-2539 FAX: (508) 845-9258 www.profoldin.com info@profoldin.com


INSTRUCTIONS

ProFoldin MicroMolar UDP Assay Kits

MicroMolar UDP assay kit Catalog number: MUD100K MicroMolar UDP assay kit - 500 Catalog number: MUD500K

INTRODUCTION

The MicroMolar UDP Assay is for measurement of UDP concentrations in a sub-micromolar to low micromolar range. Other nucleoside diphosphates (ADP, GDP and CDP) showed very little background. Nucleoside triphosphates (ATP, UTP, GTP and CTP) and monophosphates (AMP, UMP, GMP and CMP) are not detected. The UDP assay is based on fluorescence measurement with emission at 535 nm and excitation at 485 nm. It is a high throughput assay using regular black 96-well or 384-well plates. Microcuvettes may also be used for detection of the fluorescence signals. The assay can be used to measure contamination of UDP in UTP samples or monitor hydrolysis of UTP. It can also be used for measurement of enzyme activities that generate UDP molecules. The UDP assay is compatible with common reaction buffers with magnesium. It is not compatible with samples containing high concentrations of DNA or RNA.

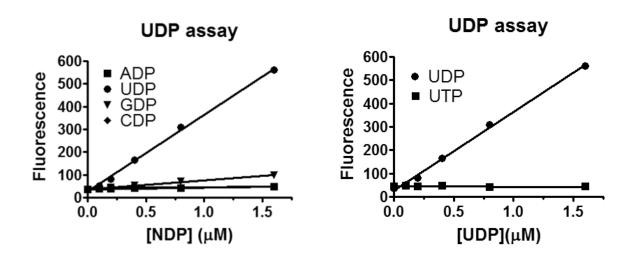

MicroMolar UDP assay kit (Catalog number: MUD100K): The kit contains 350 μl of 10 x Buffer, 30 μl of 100 x MUD reagent 1, 30 μl of 100 x MUD reagent 2, and 300 μl of 10 x fluorescence dye. The assay reagents are sufficient for measurement of 100 samples of UDP using a standard black 384-well plate.

MicroMolar UDP assay kit - 500 (Catalog number: MUD500K): The kit contains $1800 \,\mu l$ of $10 \,x$ Buffer, $150 \,\mu l$ of $100 \,x$ MUD reagent 1, $150 \,\mu l$ of $100 \,x$ MUD reagent 2 and $1500 \,\mu l$ of $10 \,x$ fluorescence dye. The assay reagents are sufficient for measurement of $500 \,samples$ of UDP using a standard black 384-well plate.

ProFoldin

10 Technology Drive, Suite 40, Number 188 Hudson, MA 01749-2791 USA Tol. (508) 735-2530 FAX: (508) 845-025

Tel: (508) 735-2539 www.profoldin.com FAX: (508) 845-9258 info@profoldin.com

INSTRUCTIONS

PROTOCOL

The following assay protocol is based on using a 384-well plate (Matrix 4318 plate) for the measurement. The sample volume is 15 μ l and the final assay volume is 30 μ l. For 96-well plate assays, the sample volume is 50 μ l and the final assay volume is 100 μ l. For assays using cuvette, the sample volume is 500 μ l and the final assay volume is 1000 μ l.

STANDARD CURVE

1. **Sample preparation:** Prepare 15 μl of UDP solutions in a 384-well black plate (Matrix 4318 plate) with a two-fold serial dilution from 0.005 mM to zero in water. Freshly prepare the following premix for each 10 assays: 136.8 μl of H₂O, 36 μl of 10 x Buffer, 3.6 μl of 100 x MUD Reagent 1, 3.6 μl of 100 x MUD Reagent 2.

The 1 x Buffer is composed of 50 mM Tris-HCl, pH 8.0, 3 mM MgCl₂, 0.2 mM EDTA, 0.5 mM DTT, 50 mM NaCl, 0.003% Brij-35.

- 2. **Detection:** Mix 15 μl of the UDP samples with 15 μl of the premix for 45 min. Dilute 33μl of the 10 x Fluorescence dye with 297 μl of water (10-fold dilution). Add 30 μl of 1 x Fluorescence dye and read the fluorescence intensity at 535 nm with excitation at 485 nm in 5 min.
- 3. **Data Analysis**: Plot the fluorescence intensity **Fc** and the UDP concentration [**UDP**] to generate the linear standard curve.

$$Fc = a [UDP] + b$$

Where the **Fc** values are from experimental data, the **a** and **b** values are from the linear fitting between the **Fc** values and the UDP concentrations.

UNKNOWN SAMPLES

Follow the same procedure to measure the fluorescence intensity \mathbf{Fc} values from the unknown samples. Calculate the UDP concentrations in the unknown samples using the \mathbf{Fc} values from the unknown samples and the \mathbf{a} and \mathbf{b} values from the standard curve.

$$[UDP] = (Fc - b) / a$$

PUBLICATIONS

- 1. Takayama F. et al, Diurnal dynamic behavior of microglia in response to infected bacteria through the UDP-P2Y6 receptor system. Sci Rep. 6: 30006 (2016).
- 2. Sobhanifar S. et al, Structure and mechanism of Staphylococcus aureus TarM, the wall teichoic acid α-glycosyltransferase, PNAS, 112 (6) E576-E585 (2015).

RELATED PRODUCTS

MicroMolar ADP assay kit

MicroMolar ADP assay kit - 500

MicroMolar GDP assay kit

MicroMolar GDP assay kit

MicroMolar GDP assay kit - 500

MicroMolar Universal Kinase Assay Kit-500 (kinase not included)

Thymidylate Kinase Assay Kit Plus-500 (kinase included)

Human Thymidylate Kinase Assay Kit Plus-500 (enzyme included)

Catalog number: MGD100K

Catalog number: MGD500K

Catalog Number: MUK500K

Catalog number: TMK500KE

Catalog number: TMK500KE