# ProFoldin 10 Technology Drive, Suite 40 Hudson, MA 01749-2791 USA Tel: (508) 735-2539 FAX: (508) 845-9258 www.profoldin.com info@profoldin.com

### INSTRUCTIONS

# **ProFoldin**

# E. coli UMP Kinase Assay Kit Plus


E. coli UMP Kinase Assay Kit Plus-100 E. coli UMP Kinase Assay Kit Plus-500 Catalog number: UMK100KE Catalog number: UMK500KE

#### Introduction

UMP kinase or uridylate kinase is essential for biosynthesis of nucleic acid. It catalyzes the ATP-dependent phosphorylation of UMP into UDP.

$$UMP + ATP \longrightarrow UDP + ADP$$

UMP kinase is an attractive antibacterial target. The UMP kinase assay is based on measurement of ADP generated from the kinase reaction. The assay is fluorescence-based and can be carried out using regular black 96-well or 384-well plates or micro-cuvettes.

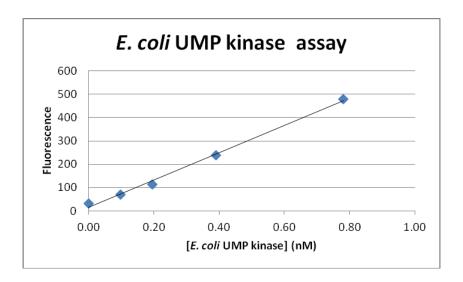

The *E. coli* UMP Kinase Assay Kit Plus (Catalog No. UMK100KE) includes 600 μl of 10 x assay buffer, 35 μl of 100 x UMP, 35 μl of 100 x *E. coli* UMP kinase, 35 μl of 100 x ATP, 35 μl of 100 x MUK Reagent A, 35 μl of 100 x MUK Reagent B and 300 μl of 10 x fluorescence dye for 100 assays in a 384-well assay format.

The *E. coli* UMP Kinase Assay Kit Plus-500 (Catalog No. UMK500KE) includes 2 ml of 10 x assay buffer, 170 μl of 100 x UMP, 170 μl of 100 x *E. coli* UMP kinase, 170 μl of 100 x ATP, 170 μl of 100 x MUK Reagent A, 170 μl of 100 x MUK Reagent B and 1.7 ml of 10 x fluorescence dye for 100 assays in a 384-well assay format.

# ProFoldin 10 Technology Drive, Suite 40 Hudson, MA 01749-2791 USA Tel: (508) 735-2539 FAX: (508) 845-9258 www.profoldin.com info@profoldin.com

## **INSTRUCTIONS**

#### **Assay Protocol**

#### 1. Reagent preparation:

- (1) 1 x reaction buffer: dilute the 10 x Buffer 10-fold with water. The 1 x buffer is composed of 50 mM Tris-HCl, pH 8.0, 3 mM MgCl<sub>2</sub>, 0.2 mM EDTA, 0.5 mM DTT, 50 mM NaCl, 0.003% Brij-35.
- (2) 10 x UMP: dilute the 100 x UMP solutions (10 mM) 10-fold with water.
- (3) 10 x ATP: dilute the 100 x ATP solutions (2 mM) 10-fold with water.
- (4) 10 x UMP kinase: dilute the 100 x UMP kinase 10-fold with 1 x buffer.
- (5) 10 x MUK Reagent A: dilute the 100 x MUK Reagent A 10-fold with water.
- (6) 10 x MUK Reagent B: dilute the 100 x MUK Reagent B 10-fold with 1 x buffer.
- (7) 1 x fluorescence dye: dilute the 10 x Fluorescence dye 10-fold with water.

#### 2. Kinase assay

The kinase reaction volume is 30 µl.

- (1) In a standard black 384-well plate (Matrix 4318), mix 12 μl of water, 3 μl of 10 x buffer, 3 μl of 10 x UMP, 3 μl of 10 x ATP, 3 μl of 10 x kinase. Incubate the reaction for 2 min.
- (2) Add 3 µl of 10 x MUK Reagent A, 3 µl of 10 x MUK Reagent B.
- (3) Incubate the reaction mixture at 37°C for 60 min.
- (4) Add 30 µl of the 1 x fluorescence dye into the 30 µl of the reaction mixture.
- (5) Measure the fluorescence intensity at 535 nm with excitation at 485 nm.

### Assay Protocol for enzyme inhibition

In order to accurately measure the  $IC_{50}$ , it is important to make sure the assay is in the linear range. The linear range can be defined by enzyme concentration-dependence and time-dependence experiments. The enzyme concentration and time in the linear range is selected for the IC50 measurement.

For the enzyme inhibition experiment, typically 50 x stock solutions with a 2-fold serial dilution in water or DMSO are prepared. In the wells of the 384-well plate,  $0.6 \mu l$  of the 50 x inhibitor is mixed with a premix composed of the buffer and enzyme for 5 min. Then the substrate is added. After incubation for 2 min,  $3 \mu l$  of 10 x MUK Reagent A and  $3 \mu l$  of 10 x MUK Reagent B are added. The total assay volume is  $30 \mu l$ . At the end of the reaction, the dye is added and the fluorescence intensity is measured.

# **Related products**

E. coli Guanylate Kinase Assay Kit Plus-500
 E. coli NAD+ Kinase Assay Kit Plus-500
 E. coli Thymidylate Kinase Assay Kit Plus-500
 Catalog number: NAK500KE
 Catalog number: TMK500KE
 Human Thymidylate Kinase Assay Kit Plus-500
 Catalog number: HTMK500KE

For more information of drug targets and enzyme assays, please visit www.profoldin.com.