


Mobicol "F" with fixed outlet plug and 2 different screw caps, without filters

Product Information Sheet # M1050

SUMMARY

shipped at RT; store at RT

For research use only

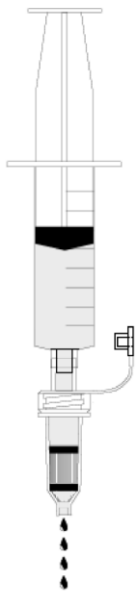

Syringe connected to Luer-lock cap


Mobicol "F" is a 1 ml column with particular handling advantages.

Mobicol "F"

- is compatible to the laboratory standard,
- can be centrifuged in minifuges,
- has a Luer-lock adapter,
- smallest and largest volumes can be treated easily.


Filters of the pore sizes 10 µm, 35 µm or 90 µm are available for Mobicols. The filters can be inserted easily with the help of a plastic tool which is delivered with the columns. The upper large filter fits tightly over the whole length of the column body.

Mobicols can be closed tightly with a screw-on cap and the reversed side of the snap-off plug at the outlet. The volume is 700 µl. Mobicols can be autoclaved at 120 °C (with inserted filter at 110 °C for 10 minutes).

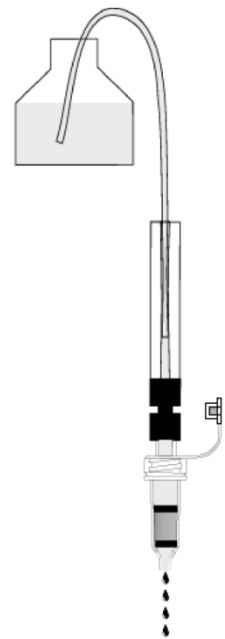

Adapter M3001: tubing to Mobicol

Mobicol "F"


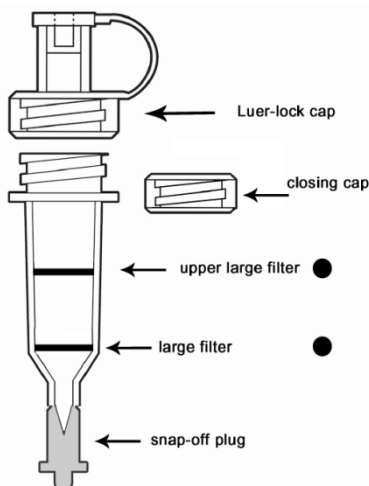


Mobicols can be centrifuged (fit into 1.5 ml or 2.2 ml tubes)

Mobicol "F" with fixed outlet plug and 2 different screw caps, without filters



Product Information Sheet # M1050

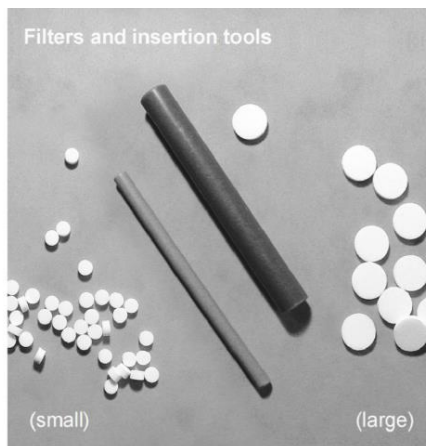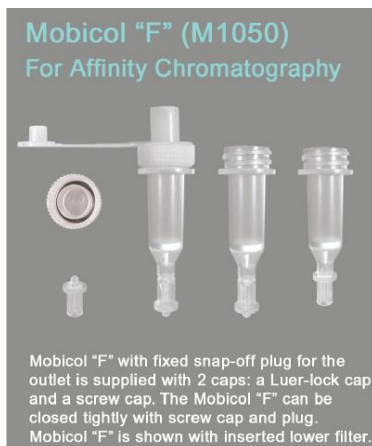

The Mobicol "F" M1050 is supplied with two screw caps (Luer-lock cap and screw cap) and a snap-off plug for the outlet. Filters with different pore sizes (10 μm , 35 μm and 90 μm) are separately available, so you can use the filter of your choice. With this features the Mobicol "F" can be used for a variety of applications.

Order Information, Shipping and Storage

| Order# | Product | Quantity |
|----------|-------------------------------------------------------------------------------------------------|----------|
| M1050 | Mobicol "F" with fixed outlet plug and 2 different screw caps, without filters | 50 |
| M1053 | Mobicol "F" with fixed outlet plug and screw cap, without filters | 50 |
| M105035F | Mobicol "F" with fixed outlet plug, inserted 35 μm filter and 2 different screw caps | 50 |
| M2210 | Filter (large) 10 μm pore size, for Mobicol M1002, M1003, M1050 & M1053 | 50 |
| M523515 | Filter (large) 35 μm pore size, for Mobicol M1002, M1003, M1050 & M1053 | 50 |
| M2290 | Filter (large) 90 μm pore size, for Mobicol M1002, M1003, M1050 & M1053 | 50 |

shipped at RT; store at RT

Column Accessories

| Order# | Product | Quantity |
|--------|----------------------------------------------------------------------------------------------------------------------------------------|----------|
| M3001 | Tubing-to-Luer-lock adapter (material outside: metal; inside: glass and PTFE; with 1 m PTFE tubing) for Mobicols or laboratory columns | 1 |
| M3002 | Luer adapter, (M+F) 10 each | 20 |
| M3003 | Luer adapter for the top-cap (M) | 20 |
| M3004 | Luer adapter for outlet (F) | 20 |
| M3006 | Stopcocks, one-way, Luer-lock | 4 |
| M3007 | Luer-lock adapter (M/M) | 1 |

shipped at RT; store at RT

Contact and Support

MoBiTec GmbH ● Lotzestr. 22a ● D-37083 Goettingen ● Germany

Customer Service – General inquiries & orders

phone: +49 (0)551 707 22 0
fax: +49 (0)551 707 22 22
e-mail: order@mobitec.com

Technical Service – Product information

phone: +49 (0)551 707 22 70
fax: +49 (0)551 707 22 77
e-mail: info@mobitec.com

MoBiTec in your area: Find your local distributor at www.mobitec.com