

Mobicol "F" with fixed outlet plug and 1 closed screw cap, without filters

Product Information Sheet # M1053

SUMMARY

shipped at RT; store at RT

For research use only

Mobicol "F" is a 1 ml column with particular handling advantages.


Mobicol "F"

- is compatible to the laboratory standard,
- can be centrifuged in microfuges,
- smallest and largest volumes can be treated easily.

Filters of the pore sizes 10 µm, 35 µm or 90 µm are available for Mobicols. The filters can be inserted easily with the help of a plastic tool which is delivered with the columns. The upper large filter fits tightly over the whole length of the column body.

Mobicols can be closed tightly with a screw-on cap and a plug at the outlet. The volume is 800 µl with a small filter and 700 µl with a large filter inserted. Mobicols can be autoclaved at 120 °C (with inserted Filter at 110 °C for 10 min).

Mobicol "F"

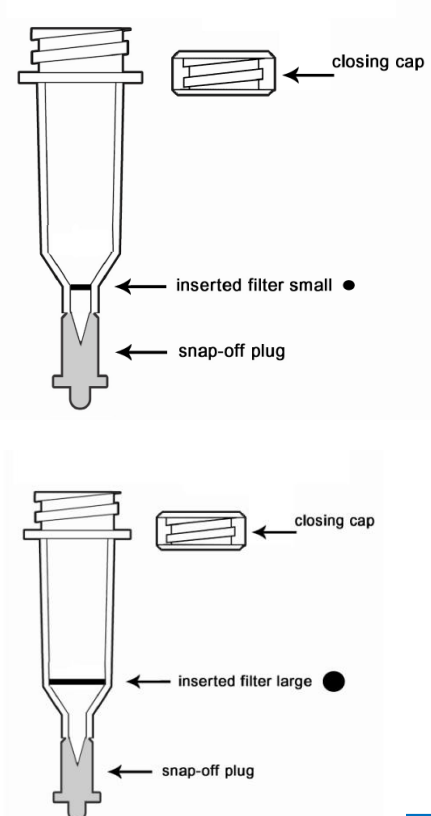


Mobicols can be centrifuged (fit into 1.5 ml or 2 ml tubes)

Mobicol "F" with fixed outlet plug and 1 closed screw cap, without filters



Product Information Sheet # M1053

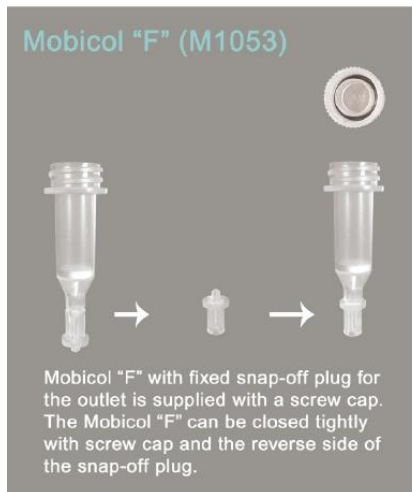

The Mobicol "F" M1053 is supplied with a screw cap and two tools for insertion of filters. The reverse side of the snap-off plug can be used for resealing. Small and large filters with different diameters (10 μm , 35 μm and 90 μm) are separately available, so you can use the filter of your choice. With this features the Mobicol "F" can be used for a variety of applications.

Order Information, Shipping and Storage

Order#	Product	Quantity
M1053	Mobicol "F" with fixed outlet plug and screw cap, without filters	50
M2110	Filter (small) 10 μm pore size, for Mobicol M1002, M1003, M1050 & M1053	50
M513515	Filter (small) 35 μm pore size, for Mobicol M1002, M1003, M1050 & M1053	50
M2190	Filter (small) 90 μm pore size, for Mobicol M1002, M1003, M1050 & M1053	50
M2210	Filter (large) 10 μm pore size, for Mobicol M1002, M1003, M1050 & M1053	50
M523515	Filter (large) 35 μm pore size, for Mobicol M1002, M1003, M1050 & M1053	50
M2290	Filter (large) 90 μm pore size, for Mobicol M1002, M1003, M1050 & M1053	50
shipped at RT; store at RT		

Contact and Support

MoBiTec GmbH ● Lotzestrasse 22a ● D-37083 Goettingen ● Germany

Customer Service – General inquiries & orders

phone: +49 (0)551 707 22 0
fax: +49 (0)551 707 22 22
e-mail: order@mobitec.com

Technical Service – Product information

phone: +49 (0)551 707 22 70
fax: +49 (0)551 707 22 77
e-mail: info@mobitec.com

MoBiTec in your area: Find your local distributor at www.mobitec.com