

Kex2 Protease (Lys/Arg-Arg), recombinant

Product Information Sheet
EP0410/EP0450



SUMMARY

shipped on dry ice; store at -80 °C

For research use only

Product Description and Application

- specific serine endoproteinase
- cleaves amino acid sequence N-Arg-Arg/-C and N-Lys-Arg/-C at the carboxyl end
- suited for protein sequencing and cleavage of fusion proteins with an appropriate recognition sequence

In the series of endoproteinases we also offer the newly developed endoproteinase Lys/Arg-Arg. The specific serine endoproteinase (universal serine-protease) has a MW of 68 kDa. It cleaves at the carboxyl end of the recognition sequences: Arg-Arg/X and Lys-Arg/X and thus provides new possibilities for e.g. the development of fusion protein systems with alternative cleavage sites.

Technical Details

Source:	<i>Saccharomyces cerevisiae</i>
Activators: pH-	Endoproteinase Lys/Arg-Arg is strongly Ca ²⁺ dependent and has a optimum at pH 7.0
Inhibitors:	Ala-Lys-Arg-chloro-methyl ketone Note: <i>Endoproteinase Lys/Arg-Arg is not inhibited by phenyl-methyl-sulfonyl-fluoride and tosyl-lysine-chloro-methylketone.</i>
Reaction buffer:	50 mM Tris-HCl or HEPES, 5 mM CaCl ₂ , pH 7.0 (0.5 mM PMSF; 0.1% Triton X-100 is not included, but might be added if required by your application). For fusion proteins the reaction conditions have to be determined empirically.
Substrate solution:	Benzoyloxycarbonyl-L-tyrosyl-L-lysyl-L-arginin-4-nitroanilid (Z-L-Tyr-Lys-Arg-pNA ° TFA salt, Bachem No. L1250). The substrate is dissolved in reaction buffer (OD ₃₁₅ = 12). This solution can be stored at -20 °C for several months. Thaw at room temperature immediately before use.
Unit definition:	1 unit endoproteinase Lys/Arg-Arg releases 1 µmole 4-nitroaniline per minute in reaction buffer at pH 7.0 at RT.
Specific activity:	2.2 U/ml

Kex2 Protease (Lys/Arg-Arg), recombinant

Product Information Sheet
EP0410/EP0450



Assays:

The substrate solution ($OD_{315} = 12$) was thawed at RT. Assay solution: 100 μ l of the substrate solution was mixed with 400 μ l reaction buffer and put into a 500 μ l cuvette (1 cm optical path length). For this solution the photometer was adjusted to 0.0 and a range from 0 to 0.5 OD_{405} was chosen. The assay solution was mixed with 10 μ l of a 1:10 dilution of the endoproteinase Lys/Arg-Arg. The optical density was measured at 405 nm for 5 minutes (at RT). The initial slope was plotted as a straight line and, after 1 minute, the value of the straight line was read (ΔOD_{405} , for 1 minute).

$$\frac{\Delta OD_{405} \times 1 \text{ min} \times 100 \times 10}{\epsilon \times 2} = \text{activity in units/min} \times \text{ml (undiluted)}$$

(Note: $\epsilon = 9.94$)

References

- Achstetter T., *Biol. Chem.*, 259 (1984) 13334-48
- Achstetter T., *EMBO*, 4, No. 1 (1985) 173-177
- Dmochowska A., *Cell*, 50 (1987) 573-584
- Fuller R. *et al.*, *Arch. of Biochem. and Biophys.*, 207 No. 2 (1981) 445-454
- Fuller R. *et al.*, *Arch. of Biochem. and Biophys.*, 226 (1983) 292-305
- Fuller R. *et al.*, *Ann. Rev. Physiol.*, 50 (1988) 345-362
- Fuller R., *Proc. Acad. Sci. USA*, 86 (1989) 1434
- Fuller R. *et al.*, *Science*, 246 (1989) 482
- Julius, *Cell*, 36 (1984) 309
- Julius, *Cell*, 37 (1984) 1075-1089
- Mizuno, *Biochem. and Biophys. Res. Commun.*, 144 No. 2 (1987) 807
- Wagner J.C., *FEBS-Letters*, 218 (1987) 31-34
- Wagner J.C., *FEBS-Letters*, 221 No. 2 (1987) 423-426

Kex2 Protease (Lys/Arg-Arg), recombinant

Product Information Sheet
EP0410/EP0450



Order Information, Shipping and Storage

Order#	Product	Quantity
EP0410	Kex2 Protease (Lys/Arg-Arg), recombinant	10 U
EP0450	Kex2 Protease (Lys/Arg-Arg), recombinant	50 U
shipped on dry ice; store at -80 °C		

Contact and Support

MoBiTec GmbH ● Lotzestrasse 22a ● D-37083 Goettingen ● Germany

Customer Service – General inquiries & orders

phone: +49 (0)551 707 22 0
fax: +49 (0)551 707 22 22
e-mail: order@mobitec.com

Technical Service – Product information

phone: +49 (0)551 707 22 70
fax: +49 (0)551 707 22 77
e-mail: info@mobitec.com

MoBiTec in your area: Find your local distributor at www.mobitec.com